

# Application Data Sheet

## No. 177

### **System Gas Chromatograph**

C6 and Lower Boiling Hydrocarbons in Olefin Free Naphthas Nexis GC-2030LBH1
GC-2014LBH1

This method is for determining the composition of olefin free Naphtha as described in below compound table. It requires the use of a dedicated gas chromatographic system which is configured with an automatic liquid injector.

#### **Analyzer Information**

#### **System Configuration:**

One SPL injector / one capillary column / one FID detector

#### Sample Information:

Determining benzene, individual hexanes and lower boiling hydrocarbons in olefin free naphthas having a final boiling point of 260°C or lower.

#### Methods met:

**UOP-551** 

#### **Concentration Range:**

No.	Name of	Concentration Range	
	Compound	Low Conc.	High Conc.
1	Propane	100 ppmwt	10,000 ppmwt
2	Isobutane	100 ppmwt	10,000 ppmwt
3	n-Butane	100 ppmwt	10,000 ppmwt
4	Isopentane	100 ppmwt	50 %wt
5	n-Pentane	100 ppmwt	40 %wt
6	2,2-Dimethylbutane	100 ppmwt	10 %wt
7	Cyclopentane	100 ppmwt	10,000 ppmwt
8	2,3-Dimethylbutane	100 ppmwt	10 %wt
9	2-Methylpentane	100 ppmwt	40 %wt
10	3-Methylpentane	100 ppmwt	30 %wt
11	n-Hexane	100 ppmwt	20 %wt
12	Methylcyclopentane	100 ppmwt	10 %wt
13	Benzene	100 ppmwt	10 %wt
14	Cyclohexane	100 ppmwt	10,000 ppmwt

Detection limits may vary depending on the sample. Please contact us for more consultation.

#### **System Features**

- ·Single FID channel
- Good repeatability

#### **Typical Chromatograms**

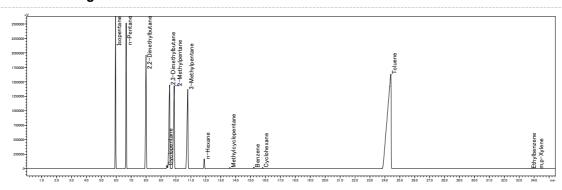


Fig. 1 Chromatogram of FID

to change without notice.

First Edition: November, 2017